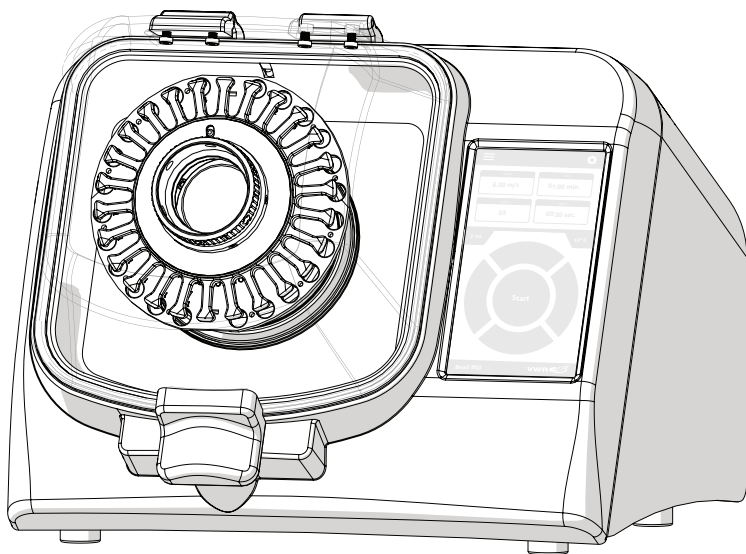
Bead Mill MAX

Bead Mill Homogeniser

User Manual

| | |
|----|----------------------|
| en | User Manual |
| fr | Manuel d'utilisation |
| it | Manuale d'uso |
| es | Manual del usuario |
| de | Benutzerhandbuch |
| po | Manual do usuário |

| Catalog No | Description |
|------------|-----------------------|
| 432-0380 | Bead Mill, 230V, 300W |

Legal Address of Manufacturer

| United States | Europe |
|---|---|
| VWR International, LLC | VWR International bv |
| 100 Matsonford Rd | Researchpark Haasrode 2020 |
| Radnor, PA 19087 | Geldenaaksebaan 464 |
| 800-932-5000 | 3001 Leuven |
| http://www.vwr.com | + 32 16 385011 |
| | http://be.vwr.com |

Country of Origin

Made in USA


Intended Use

The VWR Bead Mill is intended for milling and homogenising of a number of substances. Not intended for medical or diagnostics purposes.

Table of Contents

- 4** Overview / Package Contents
Voltage
- 5** Specifications
Loading Sample Tubes
- 6** General Operation
Overview
Run Time
- 7** Settings Menu
Language Setting
- 8** Programming Mode
- 9** Troubleshooting
- 10** Technical Service
Warranty
Equipment Disposal
- 11** Local VWR Offices in Europe and Asia Pacific

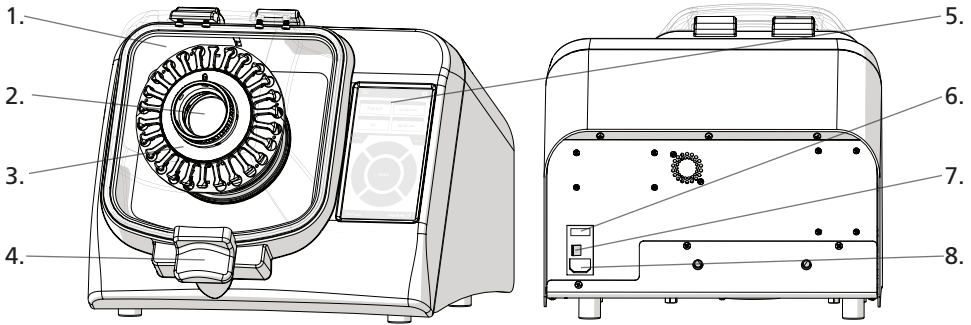Overview / Package Contents


- | | |
|------------------|----------------------------|
| 1. Lid | 5. Touch Screen |
| 2. Drive Hub | 6. 120V/220V Switch & Fuse |
| 3. Tube Carriage | 7. ON/OFF Switch |
| 4. Lid Latch | 8. Power Cord Port |

The VWR Bead Mill consists of the following:

| Description | Quantity |
|--------------------|----------|
| Bead Mill Assembly | 1 |
| Power Cord | 1 |
| Spare Fuse | 1 |
| User Manual | 1 |

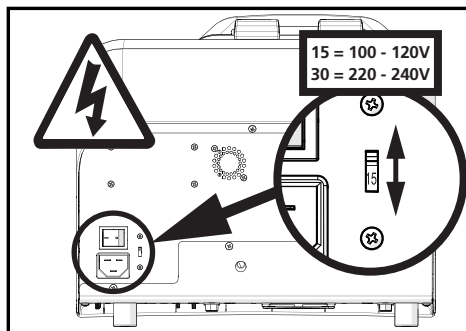Voltage

Remove Warning stickers from the back of the unit and check that the voltage matches the voltage delivered by your electric power supply.



CAUTION: DO NOT connect the unit to the main power supply until correct voltage is selected. Plug the unit into power supply using the supplied power cord.

CAUTION: If the voltage supplied to the unit is higher than what is indicated on the voltage switch, the electronics can be irreversibly damaged, voiding the warranty.



Specifications

| | |
|-------------------|--|
| Processing Range | 25 micro-liters to 50 mL |
| Power Rating | 100-230 VAC/ 50-60Hz |
| Dimensions | L: 15in(38.1cm), W: 17in(43.18cm), H: 13in(33.0cm) |
| Product Weight | 65 lbs. (29kg) |
| Compliance | EU CE Legislation |
| Run Time | 1 second to 9:59 minutes |
| Performance Range | 0.8m/s – 6.5m/s in increments of 0.15m/s |
| Number of Cycles | 1-10 |

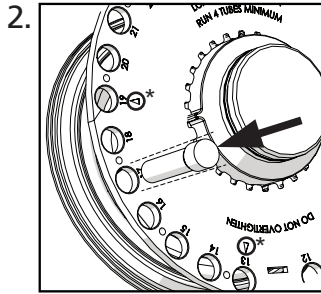
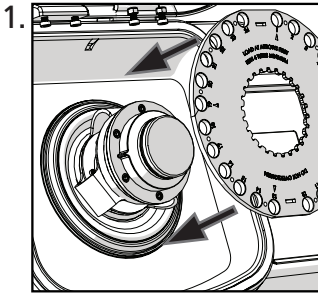
Loading Sample Tubes

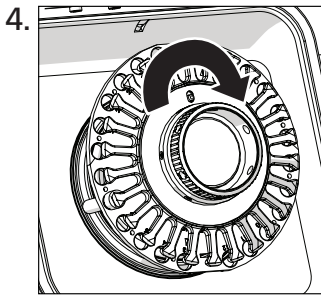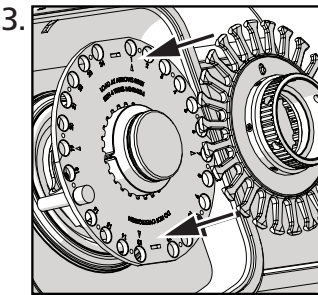A minimum of four tubes are required and must be spaced evenly*. (See figure 2)

Screw down to secure tubes.

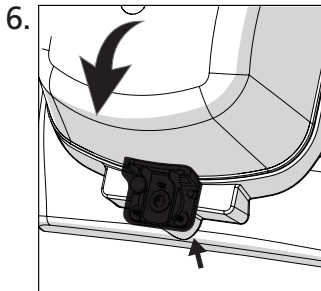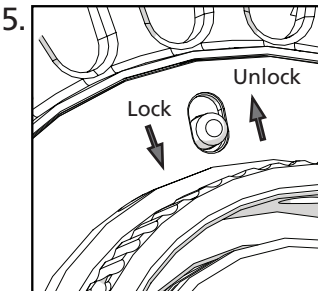Move the locking switch to the "LOCK" position.

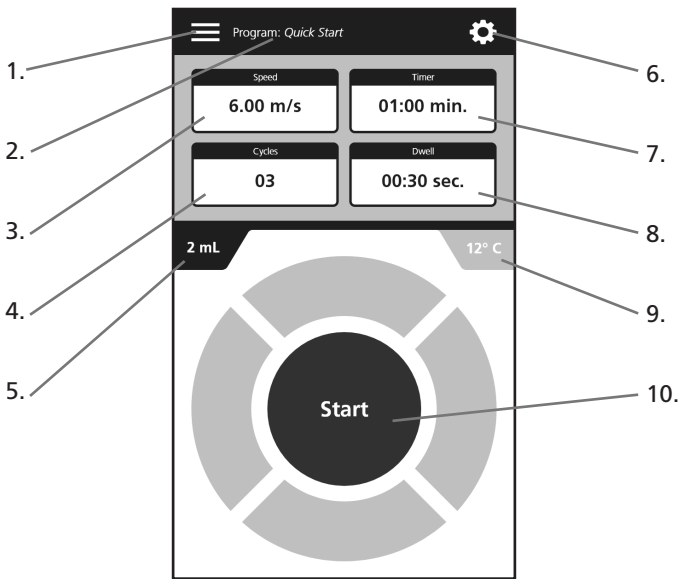
CAUTION: DO NOT run without moving the switch to the "Lock" position.

Close the lid until the locking handle clicks into place.

CAUTION: The unit will not run without the lid fully closed.


General Operation

Overview


- | | |
|---------------------|---|
| 1. Program Menu | 6. Settings Menu |
| 2. Program Name | 7. Timer |
| 3. Speed | 8. Dwell / Time Interval between Cycles |
| 4. Number of Cycles | 9. Processing Chamber Temperature |
| 5. Tube Setting | 10. Start / Select / Pause Button |

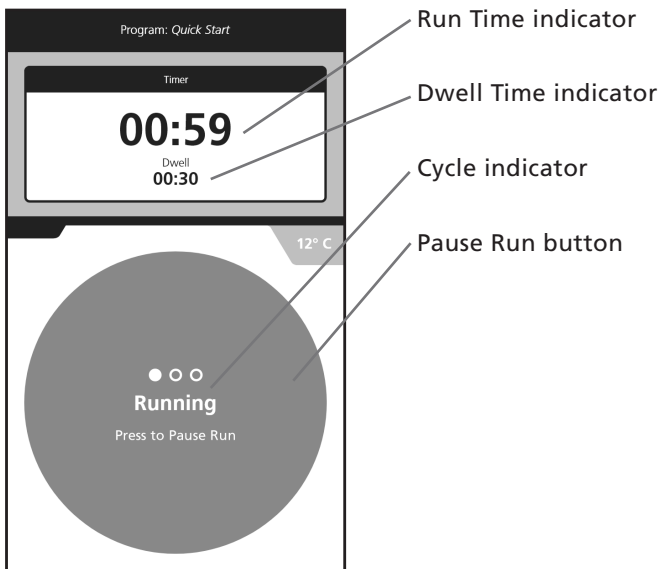Run Time


General Operation

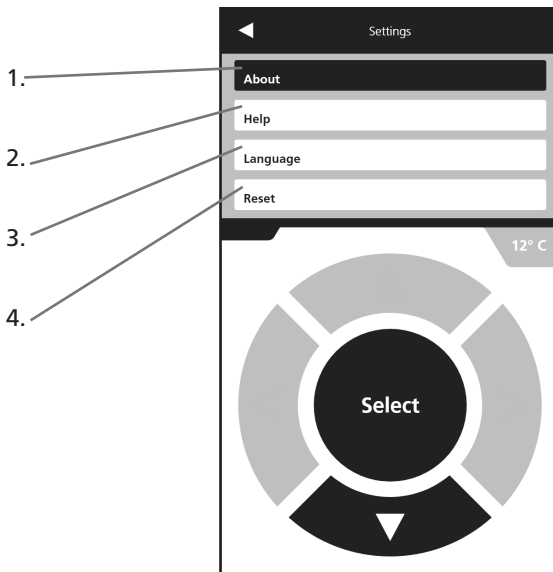Settings Menu

1. About
2. Help
3. Language
4. Reset

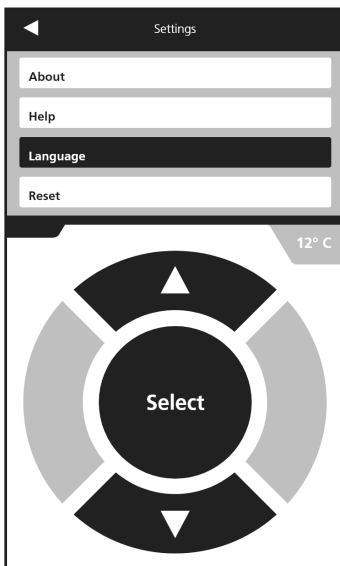
Language Setting

- Select Settings (⚙️)
- Select "Language"
- △▽ to navigate through menu
- Choose from French, English, German, Italian, or Spanish
- Press "Select"

General Operation

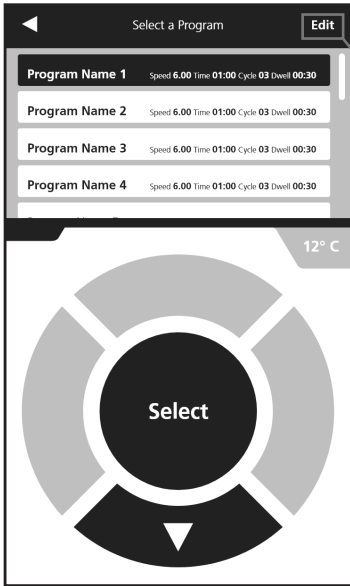Programming Mode

- Select Program Menu (☰)

△▽ to select a program from list

Press "Edit" to edit selected program

- Select variable with △▽

- Press "Edit" to adjust variable

◀▶ to select values for Time, Cycle, and Dwell

Press "Save" to save your program

| | Parameter | Operating Range |
|---|-------------------------------|---|
| S | Speed / Energy | From 0.8 m/s to 6.50 m/s |
| T | Time | From 0:01 to 9:59 in increments of 1 second |
| C | Number of Cycles | 1 to 10 |
| D | Dwell (Pause between Runs) | 0:01 to 9:59 |

Troubleshooting

| Error | Cause | Possible Solution |
|---|---|--|
| Unclosed Lid | The lid is not latched properly | 1. Check that nothing prevents the lid from closing. 2. Press lid and ensure the handle is locked properly. |
| | Detection system is faulty | 1. Turn off the unit. 2. Contact technical assistance. |
| Fan not working | No Power | 1. Check main voltage 2. Check that the voltage of the unit matches voltage being delivered by the main power supply. 3. Make sure the unit is plugged in properly. 4. Check and replace fuses if blown. |
| | Faulty Fan | 1. Turn off the unit. 2. Contact technical support. |
| Display screen shows incomplete or unusual characters | CPU needs to be reset | 1. Turn off the unit. 2. Wait 10 seconds. 3. Turn the unit back on. |
| No display on the screen | No Power | 1. Check main voltage. 2. Check that the voltage of the unit matches voltage being delivered by the main power supply. 3. Make sure the unit is plugged in properly. 4. Check and replace fuses if blown. |
| | Faulty display screen | 1. Turn off the unit. 2. Contact technical support. |
| One or several tubes are not sealed tightly | The cap is not properly sealed or the tube is faulty | If dangerous or potentially dangerous sample is contained in the tube, apply the proper decontamination procedure. |
| Unit powers on with a loud "BANG" | Motor control device damaged due to incorrect voltage selection | Return unit to VWR for replacement of motor control device. |
| Unit powers on but motor does not turn | Incorrect voltage selected | Power off the unit and select appropriate voltage. If problem persists, contact VWR for assistance. |
| Fingerplate lock is difficult to disengage. | Knob position | Turn knob slightly counter-clockwise to release tension. |
| Tube leakage | Speed is too high, or high temperature or pressure inside tubes | Reduce processing time or speed |

WARNING: DO NOT attempt to service the VWR Bead Mill in a manner other than those discussed in this manual. For any issue that is unsuccessfully corrected using this guide, please contact your authorized dealer or call your local VWR office.

Technical Service

Web Resources

Visit the VWR website at www.vwr.com for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact Us

For information or technical assistance, contact your local VWR representative or visit www.vwr.com.

Warranty

VWR warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with Local Laws and Regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it is your responsibility to correctly dispose of your equipment at life cycle end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.


Thank you.

Accessories and Consumables
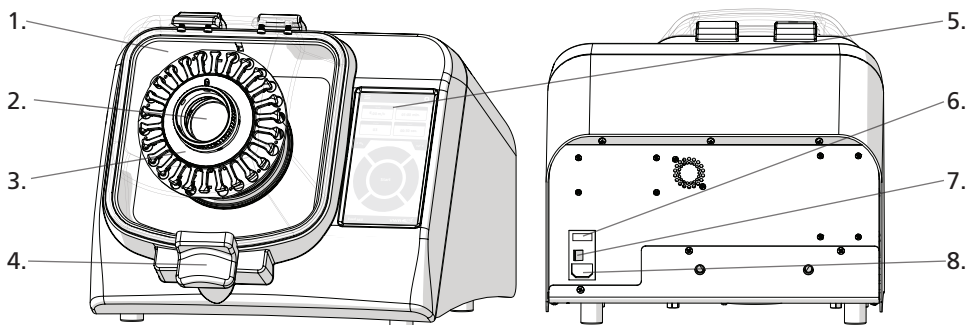
| EU Catalog No | US Catalog No | Description |
|---------------|---------------|--|
| 432-0368 | 10158-598 | 1.5 ML SCREW CAP REINFORCED TUBES PRE-FILLED 2.4 MM METAL BEADS. 50 PER PACK. |
| 432-0369 | 10158-600 | 1.5 ML SCREW CAP TUBES PRE-FILLED 1.4 MM CERAMIC BEADS. 50 PER PACK. |
| 432-0370 | 10158-602 | 1.5 ML SCREW CAP REINFORCED TUBES PRE-FILLED 2.8 MM CERAMIC BEADS. 50 PER PACK. |
| 432-0371 | 10158-606 | 2 ML SCREW CAP TUBES PRE-FILLED 0.5 MM GLASS BEADS. 50 PER PACK. |
| 432-0372 | 10158-610 | 2 ML SCREW CAP TUBES PRE-FILLED 1.4 MM CERAMIC BEADS. 50 PER PACK. |
| 432-0373 | 10158-612 | 2 ML SCREW CAP REINFORCED TUBES PRE-FILLED 2.8 MM CERAMIC BEADS. 50 PER PACK. |
| 432-0374 | 10158-604 | 2 ML SCREW CAP REINFORCED TUBES PRE-FILLED 2.4 MM METAL BEADS. 50 PER PACK. |
| 432-0375 | 10158-608 | 0.5 ML SCREW CAP TUBES PRE-FILLED 1.4 MM CERAMIC BEADS. 50 PER PACK. |
| 432-0376 | 10158-550 | 2.4 MM STAINLESS STEEL BEADS BULK PACKAGED, 500G. |
| 432-0377 | 10032-374 | 1.4 MM CERAMIC BEADS BULK PACKAGED, 325G |
| 432-0378 | 10032-372 | 2.8 MM CERAMIC BEADS BULK PACKAGED, 325G |
| 432-0379 | 10158-556 | 2 ML TUBES WITH SCREW CAPS AND O-RINGS BULK PACKAGED. 1000 PER PACK. |
| 432-0394 | 19-8005 | CRYO COOLING UNIT, COMPATIBLE WITH THE BEAD MILL MAX |
| 432-0395 | 19-345-007 | 7ML TUBE CARRIAGE KIT, COMPATIBLE WITH THE BEAD MILL MAX |
| 432-0396 | 10032-384 | 30ML TUBE CARRIAGE KIT, COMPATIBLE WITH THE BEAD MILL MAX |
| 432-0397 | 10032-368 | 7ML HARD TISSUE GRINDING MIX, 50 PACK |
| 432-0398 | 10032-366 | 7ML SOFT TISSUE HOMOGENIZING MIX, 50 PACK |
| 432-0399 | 10032-364 | 7ML HARD TISSUE HOMOGENIZING MIX, 50 PACK |
| 432-0400 | | 7ML SCREW CAP TUBES, 500 PER PACK |
| 432-0401 | | 7ML SCREW CAP TUBES, 1000 PER PACK |
| 432-0402 | 10032-522 | 30ML HARD TISSUE GRINDING MIX, 50 PACK |
| 432-0403 | 76000-782 | 30ML SOFT TISSUE HOMOGENIZING MIX, 50 PACK |
| 432-0404 | 10015-242 | 30ML HARD TISSUE HOMOGENIZING MIX, 50 PACK |
| 432-0405 | | 30ML SCREW CAP TUBES, 100 PER PACK |
| 432-0406 | | 30ML SCREW CAP TUBES, 500 PER PACK |

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VWR International AB
VWR International AG
VWR International Ltd